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The left photograph shows the 'Mode' screen of the Pro-face HMI. It features a 'Menu' button in the top left, a 'Page' button with left and right arrow icons in the top center, and 'Back' and 'Next' buttons in the top right. The main title is 'Mode'. Below it, the text 'Control Mode — User Level' is displayed. The screen is divided into two columns of buttons: the left column has 'Auto' (yellow), 'Manual' (red), and 'Int. Cancel' (red); the right column has 'Operator' (yellow), 'Maintain' (yellow), 'Admin.' (red), and a 'Password' field with a '0' digit. The bottom of the screen has a dark background.

The right photograph shows the 'Flow' screen of the Pro-face HMI. It features a 'Menu' button in the top left, a 'Page' button with left and right arrow icons in the top center, and 'Back' and 'Next' buttons in the top right. The main title is 'Flow'. Below it, the text 'N₂' is displayed. The screen shows a process diagram with various components and their settings. On the left, there are two sections: 'Ar' with 'Set' (2.8sccm) and 'Mon.' (0.3sccm) values, and 'Oz' with 'Set' (0.0sccm) and 'Mon.' (2.6sccm) values. In the center, there are two 'RFG' (Refrigerant Flow Gauge) components. On the right, there are two 'PG' (Pressure Gauge) components. At the bottom, there are 'Process Timer' settings for 'FWD' and 'REF' (both 0W), and a 'Set' button. The bottom of the screen has a dark background.

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